


KKURABO

DA-c-7

Automated Genomic DNA Extraction from Nail

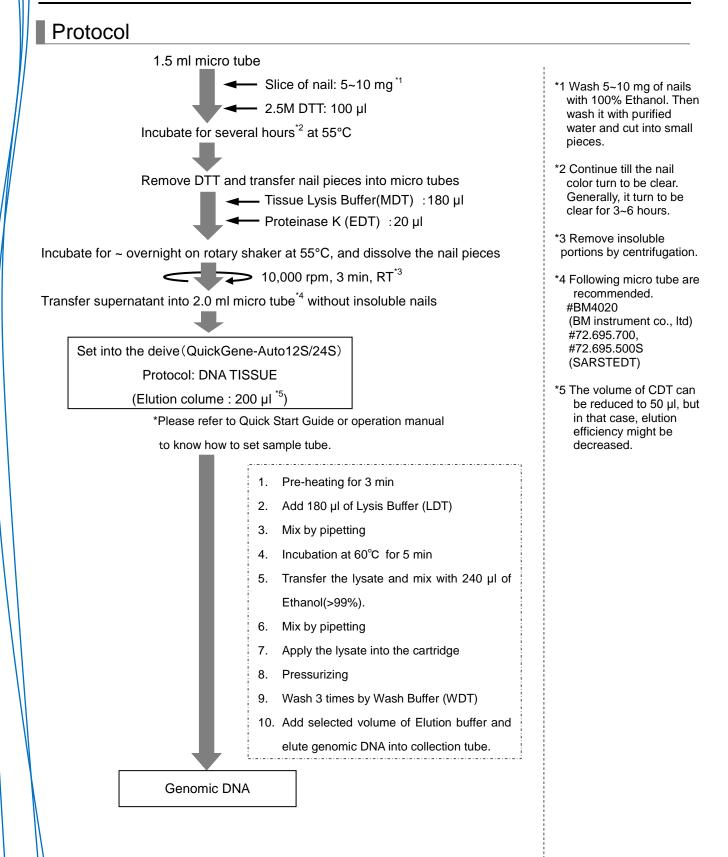

Depending on sample and storage conditions, nucleic acid may not be extractable. Therefore, we cannot guarantee accurate data. The extracted nucleic acid contains unintended acid (ex: when extracting DNA, RNA is also extracted).



Results

The yield of genomic DNA

| Sample ID | #1 | #2 | #3 | #4 | #5 | #6 |
|-------------------------------|----------|---------|---------|---------|---------|---------|
| Retention period of sample | 10 years | 9 years | 8 years | 9 years | 3 years | 3 years |
| Yield (µg) | 0.34 | 1.98 | 1.86 | 0.44 | 2.74 | 3.26 |

Protein contamination: A260/280

| SampleID | #1 | #2 | #3 | #4 | #5 | #6 |
|----------|------|------|------|------|------|------|
| A260/280 | 1.54 | 1.83 | 1.63 | 1.72 | 1.94 | 1.79 |

Common protocol is usable for the following

Depending on sample and storage conditions, nucleic acid may not be extractable. Therefore, we cannot guarantee accurate data. The extracted nucleic acid contains unintended acid (ex: when extracting DNA, RNA is also extracted).

