


DA-a-6

Automated_Genomic DNA Extraction from Whole Blood of Human

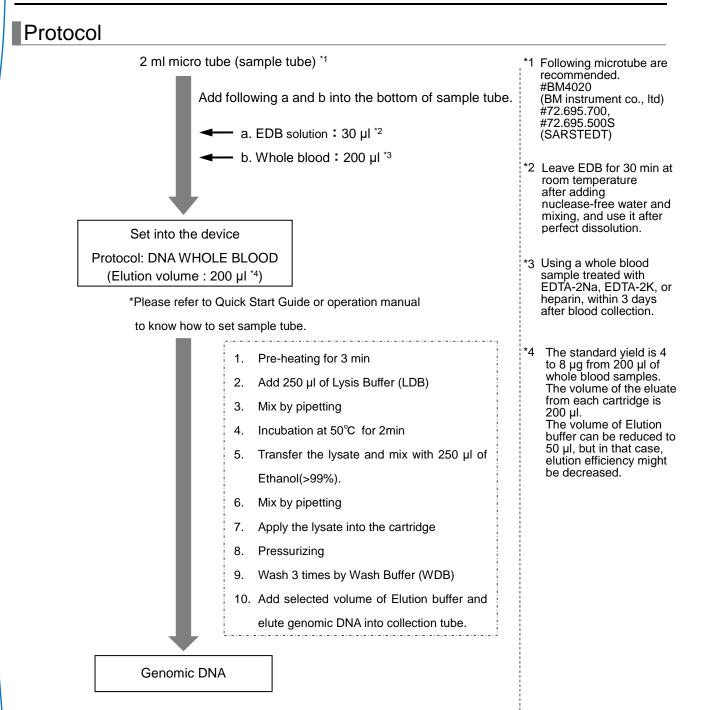

KKURABO

Depending on sample and storage conditions, nucleic acid may not be extractable. Therefore, we cannot guarantee accurate data. The extracted nucleic acid contains unintended acid (ex: when extracting DNA, RNA is also extracted).



Results

The yield of genomic DNA

	5					
Sample ID	#1	#2	#3	#4	Average	S.D.
Yield (µg)	5.6	5.4	5.6	5.5	5.5	0.10

Protein contamination : A260/280

Sample ID	#1	#2	#3	#4	Average
A260/280	1.94	2.13	2.04	2.10	2.05

Chaotropic salt contamination : A260/230

Sample ID	#1	#2	#3	#4	Average
A260/230	1.79	2.07	1.89	1.99	1.93

Common protocol is usable for the following

Canine Whole Blood

